



PEmax Enzyme

Manual

Catalog #	3473	3476
Package Size	80 µg	400 µg



Important!

-20°C Storage Required

- * Immediately inspect packages
- * Freeze upon receipt



visit us online for more
products & custom services

Intact Genomics, Inc.

Table of Contents

Product Description.....	3
Product Source.....	4
Components and Storage.....	4
Reaction Buffer.....	4
Quality Control.....	4
Functional Testing.....	5
Related Products.....	6
Ordering Information.....	6
References.....	6
Technical Support.....	7

Description:

CRISPR-associated (Cas) systems

Clustered regularly interspaced short palindromic repeats (CRISPR)/CRISPR-associated (Cas) systems provide bacteria and archaea with adaptive immunity against viruses and plasmids by using CRISPR RNAs (crRNAs) to guide the silencing of invading nucleic acids¹. The CRISPR system consists of a short noncoding guide RNA (sgRNA) made up of a target complementary CRISPR RNA (crRNA) and an auxiliary transactivating crRNA (tracrRNA)². Prime-editing is gaining popularity as a precision gene-editing technique. Prime-editing enzymes are a fusion of the Cas9 nickase mutant (H840A) with a modified Murine Moloney Leukemia Virus-Reverse Transcriptase (MMLV-RT). One of the resulting optimized prime-editing fusions is known as PEmax enzyme³⁻⁴. This prime-editing enzyme has enabled precise gene-editing without the need for double-stranded DNA breaks (DSBs) or donor DNA templates. PEmax requires a unique long gRNA known as “pegRNA” shown schematically in Figure 1.

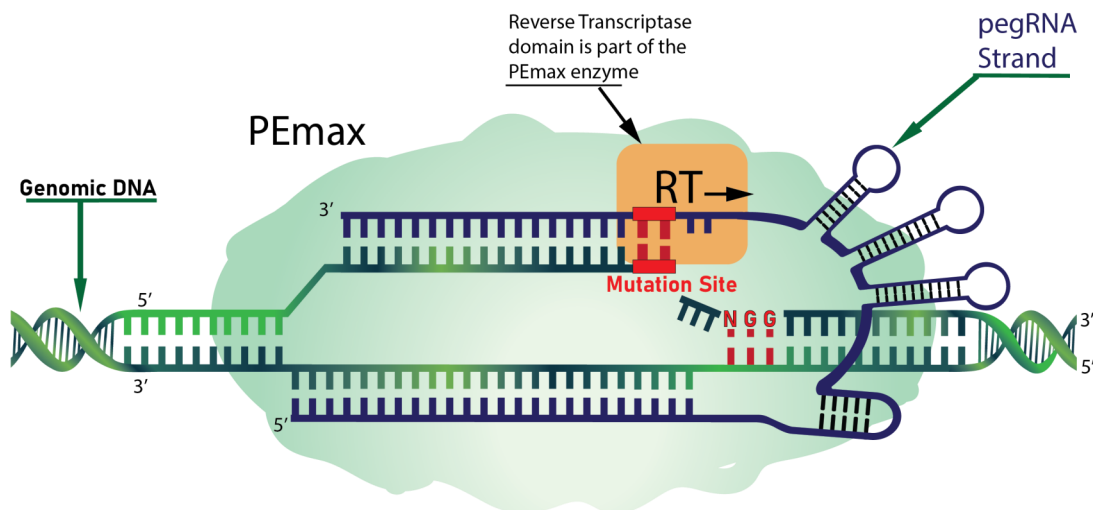


Fig. 1: Overview of the CRISPR-associated (Cas) system showing a representative prime-editing enzyme oriented with target genomic DNA and pegRNA.

Intact Genomics (IG[®]) PEmax enzyme is the purified recombinant *Streptococcus pyogenes* Cas9 nickase mutant (H840A)-Murine Moloney Leukemia Virus-Reverse Transcriptase (MMLV-RT) fusion protein³⁻⁴ containing a nuclear localization signal (NLS) at the C-terminal for targeting to the nucleus. This enzyme is designed to perform CRISPR prime-editing³⁻⁴. The physical purity of this enzyme is $\geq 85\%$ as assessed by SDS-PAGE with Coomassie[®] blue staining.

Product Source:

E. coli BL21 (DE3) strain expressing a recombinant PEmax gene from *Streptococcus pyogenes* with custom protein purification tags.

Components and Storage:

PEmax Enzyme Kit contains the items below. Store all components at -20°C.

- PEmax enzyme
- 10x PEmax Reaction Buffer

Storage Buffer:

- 50 mM Tris-HCl, 50 mM KCl, 1 mM DTT, 0.1 mM EDTA, 50% Glycerol, pH 7.5 @ 25 °C

1x Cas9 Reaction Buffer:

- 20 mM HEPES, 100 mM NaCl, 5 mM MgCl₂, 0.1 mM EDTA, pH 6.5 @ 25 °C

Quality Control:

PEmax enzyme is free from detectable RNase, Endonuclease (nicking) and non-specific DNase activities.

Functional Testing:

We have demonstrated that PEmax has specific nicking activity in the presence of an *in vitro* transcription-synthesized pegRNA guide molecule (Fig.2) and *in vitro* RT activity (not shown).

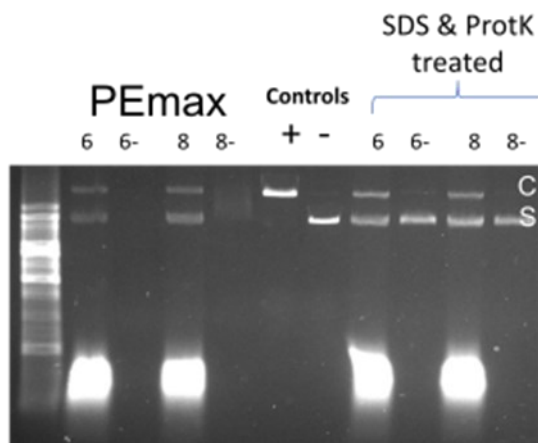


Figure 2. PEmax successfully nicks DNA, and it requires pegRNA for function. The PEmax nicking assay shows circular (C) and supercoiled (S) dsDNA. The positive (+) control uses Nb.BssSI, a nicking restriction enzyme from NEB, which nicks DNA so that it runs mostly as circular DNA (lane +). The negative (–) control is dsDNA with NO enzyme added. The DNA is mostly supercoiled; the circular DNA is too faint to see (lane –). PEmax is unable to nick without pegRNA (lanes 6- and 8-). However, in the presence of pegRNA, PEmax nicks dsDNA generating more circular dsDNA (lanes 6 and 8). For better visibility of resulting nicked DNA, SDS and Proteinase K were added to the reaction sample to remove proteins so that more DNA could enter the gel.

Related Products:

- IG® PEmax pegRNA synthesis kit (Cat.# 3403, 3406)
- IG® RNA Cleanup Kit (Cat.# 4003, 4005)
- IG® sgRNA Synthesis Kit for Cas9 Nuclease (Cat.# 3203, 3206)
- IG® sgRNA Synthesis Kit for Cas12a Nuclease (Cat.# 3303, 3306)
- Cas9 Nuclease (Cat.#3273, 3276)
- Cas12a Nuclease (Cat.# 3373, 3376)
- ig® 10B Chemically Competent Cells (Cat.# 1011-12)

Ordering Information:

- Order online within the USA. Place orders on **www.intactgenomics.com** using our secure Shopping Cart.
- Order by email, phone, or fax.
Email: **sales@intactgenomics.com**
Phone: (314) 942-3655 | Toll-free : 855-835-7172 | Fax: (314) 942-3656
- Order via our distributors.

References:

1. Jinek, M. et al. A programmable dual RNA-guided DNA endonuclease in adaptive bacterial immunity. *Science* 337, 816–821 (2012).
2. Mali, P. et al. RNA-Guided Human Genome Engineering via Cas9. *Science* 339, 823–826 (2013).
3. Chen, P. J. et al. Enhanced prime editing systems by manipulating cellular determinants of editing outcomes. *Cell* 184, 5635-5652.e29 (2021).
4. Anzalone, A. V. et al. Search-and-replace genome editing without double-strand breaks or donor DNA. *Nature* 576, 149–157 (2019).

Intact Genomics owns the following registered trademarks granted by the United States Patent and Trademark Office (USPTO): Intact Genomics®, IG®, ig®, igTherapeutics®, FastAmp®, i7®, DirectPlate®.

All technology protocols discussed within this manual are assumed proprietary to Intact Genomics. This Product may be covered by pending or issued patents or may have certain limitations. Please contact us for more information. Purchase of this material conveys to buyer the non-transferable right to use the material purchased in research conducted by buyer, whether for teaching, non-commercial or commercial research purposes. Buyer may not sell or otherwise transfer these materials, its components, or unmodified descendants to a third party.

Product Use Limitation and Disclaimers

This product is for research purposes only. It is not intended for therapeutic or diagnostic purposes in humans or animals. This product contains chemicals which may be harmful if misused or direct human contact is made.

Intact Genomics is dedicated to practicing and maintaining science and technology ethics. Buyer agrees to use the purchased materials in full compliance with applicable law and regulations.

Technical Support & Customer Services

Intact Genomics (IG®) is dedicated to customer satisfaction regarding the use of our products for your research needs. Each new lot of our products is thoroughly tested to ensure it meets high quality standards and provides excellent results. We appreciate your business and your feedback regarding the performance of our products in your applications. Please follow the instructions carefully and contact us if additional assistance is needed.

Our hours are Monday - Friday, 8AM to 5PM, U.S. Central Standard Time.

Intact Genomics, Inc.

11840 Westline Industrial Drive, Suite 120,
St. Louis, MO. 63146, USA

Phone: (314) 942-3655 | **Toll-free :** 855-835-7172 | **Fax:** (314) 942-3656

Email: sales@intactgenomics.com | ig@intactgenomics.com

Website: www.intactgenomics.com

© 2024 Intact Genomics, Inc
All Rights Reserved

