



## T4 Polynucleotide Kinase (PNK)

### Manual

<b>Catalog #</b>	<b>3232</b>
<b>Package Size</b>	2,500 units
<b>Volume</b>	250 µl
<b>Concentration</b>	10 units/µl



### Important!

#### **-20°C Storage Required**

- \* Immediately inspect packages
- \* Freeze upon receipt



visit us online for more  
products & custom services

**Intact Genomics, Inc.**

## Table of Contents

<b>Product Description</b> .....	3
<b>Physical Purity</b> .....	3
<b>Product Source</b> .....	3
<b>Applications</b> .....	3
<b>Product Components</b> .....	3
<b>Storage</b> .....	3,4
<b>Reaction Buffer</b> .....	4
<b>Unit Definition</b> .....	4
<b>Inhibition and Inactivation</b> .....	4
<b>Quality Control</b> .....	4
<b>Radioactive/Non-Radioactive Information</b> .....	4
<b>Related Products</b> .....	5
<b>Ordering Information</b> .....	5
<b>References</b> .....	5
<b>Technical Support</b> .....	6

### Description:

T4 Polynucleotide Kinase (PNK) catalyzes the transfer of the  $\gamma$ -phosphate from ATP to the 5'-OH group of single- and double-stranded DNAs and RNAs, oligonucleotides or nucleoside 3'-monophosphates. In the presence of ADP, T4 PNK exhibits 5'-phosphatase activity and catalyzes the exchange of phosphate group between 5'-P-oligonucleotides/polynucleotides and ATP (1). The enzyme is also a 3'-phosphatase (2).

### Physical Purity:

The purity of this enzyme is >98% homogeneity as determined by SDS-PAGE using Coomassie® Blue staining (see figure below).

### Product Source:

*E. coli* cells with a cloned pseT gene of bacteriophage T4.

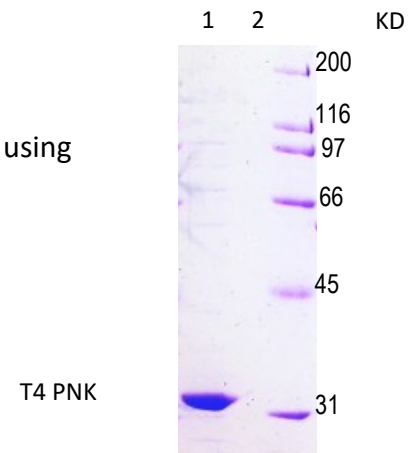


Figure: Lane 1. T4 PNK  
Lane 2. Protein Marker

### Applications:

- Labeling 5'-termini of nucleic acids (3, 4)
- 5'-phosphorylation of oligonucleotides, PCR products, other DNA or RNA prior to ligation.
- Phosphorylation of PCR primers.
- Detection of DNA modification by the [32P]-post labeling assay (5, 6).
- Removal of 3'-phosphate groups (2).

### Product Components:

- T4 Polynucleotide Kinase (PNK)
- 10x T4 PNK Reaction Buffer

### Storage Temperature:

-20 °C

### Storage Buffer:

50 mM Tris-HCl (pH 7.5), 0.1 mM EDTA, 1 mM  $\beta$ -mercaptoethanol, 1 mM DTT, 25% (v/v) glycerol

### 10x T4 PNK Reaction Buffer:

500 mM Tris-HCl, 100 mM  $MgCl_2$ , 50 mM dithiothreitol, pH 7.5 @ 25°C

### Unit Definition:

One unit of T4 Polynucleotide Kinase converts 1 nmol of  $^{32}P$  from  $[\gamma\text{-}^{32}P]\text{-ATP}$  into an acid-insoluble form in 30 minutes at 37 °C under standard assay conditions.

### Inhibition and Inactivation:

- Inhibitors: metal chelators, phosphate and ammonium ions, KCl and NaCl at a concentration higher than 50 mM.
- Inactivated by heating at 70 °C for 15 min or by addition of EDTA.

### Quality Control Assays:

T4 PNK is tested in 5' phosphorylation of nucleic acids and is free from exo- and endonuclease and RNase activities.

**Note:** T4 PNK requires ATP for activity, but it is not supplied with ATP because it interferes with radiolabeling reactions.

### For Radioactive Labeling:

Use 1–50 pmol of 5' termini in a 50  $\mu$ l reaction containing 1X T4 PNK reaction buffer, 50 pmol of gamma- $^{32}P$  ATP and 20 units of T4 PNK. Incubate at 37 °C for 60 minutes.

### For Non-Radioactive Phosphorylation:

Use up to 300 pmol of 5' termini in a 50  $\mu$ l reaction containing 1X T4 PNK reaction buffer, 1 mM ATP and 10 units of T4 PNK. Incubate at 37 °C for 60 minutes.

### Related Products:

- *Taq* DNA Polymerase (Cat.# 3243)
- *Taq* DNA Polymerase 2x Premix (Cat.# 3249)
- *Taq* DNA Ligase (Cat.# 3219)
- T4 DNA Ligase (Cat.# 3212)
- T4 DNA Polymerase (Cat.# 3222)

### Ordering Information:

- Order online within the USA. Place orders on **www.intactgenomics.com** using our secure Shopping Cart.
- Order by email, phone, or fax.  
Email: **sales@intactgenomics.com**  
Phone: (314) 942-3655 | Toll-free : 855-835-7172 | Fax: (314) 942-3656
- Order via our distributors.

### References:

1. Berkner, K.L., Folk, W.R., Polynucleotide kinase exchange reaction, *J. Biol. Chem.*, 252, 3176-3184, 1977.
2. Richardson, C.C., Bacteriophage T4 polynucleotide kinase, *The Enzymes* (Boyer, P.D., ed.), 14, 299-314, Academic Press, San Diego, 1981.
3. Sambrook, J., Russell, D.W., *Molecular Cloning: A Laboratory Manual*, the third edition, Cold Spring Harbor Laboratory Press, Cold Spring Harbor, New York, 2001.
4. *Current Protocols in Molecular Biology*, vol. 1 (Ausubel, F.M., et al., ed.), John Wiley & Sons, Inc., Brooklyn, New York, 3.10.2-3.10.5, 1994-2004.
5. Phillips, D.H., Detection of DNA modifications by the 32P-postlabelling assay, *Mutation Res.*, 378, 1-12, 1997.
6. Keith, G., Dirheimer, G., Postlabeling: a sensitive method for studying DNA adducts and their role in carcinogenesis, *Curr. Opin. Biotechnol.* 6, 3-11, 1995.

Intact Genomics owns the following registered trademarks granted by the United States Patent and Trademark Office (USPTO): Intact Genomics®, IG®, ig®, igTherapeutics®, FastAmp®, i7®, DirectPlate®.

All technology protocols discussed within this manual are assumed proprietary to Intact Genomics. This Product may be covered by pending or issued patents or may have certain limitations. Please contact us for more information. Purchase of this material conveys to buyer the non-transferable right to use the material purchased in research conducted by buyer, whether for teaching, non-commercial or commercial research purposes. Buyer may not sell or otherwise transfer these materials, its components, or unmodified descendants to a third party.

## Product Use Limitation and Disclaimers

This product is for research purposes only. It is not intended for therapeutic or diagnostic purposes in humans or animals. This product contains chemicals which may be harmful if misused or direct human contact is made.

Intact Genomics is dedicated to practicing and maintaining science and technology ethics. Buyer agrees to use the purchased materials in full compliance with applicable law and regulations.

## Technical Support & Customer Services

Intact Genomics (IG®) is dedicated to customer satisfaction regarding the use of our products for your research needs. Each new lot of our products is thoroughly tested to ensure it meets high quality standards and provides excellent results. We appreciate your business and your feedback regarding the performance of our products in your applications. Please follow the instructions carefully and contact us if additional assistance is needed.

Our hours are Monday - Friday, 8AM to 5PM, U.S. Central Standard Time.

### Intact Genomics, Inc.

11840 Westline Industrial Drive, Suite 120,  
St. Louis, MO. 63146, USA

**Phone:** (314) 942-3655 | **Toll-free :** 855-835-7172 | **Fax:** (314) 942-3656

**Email:** [sales@intactgenomics.com](mailto:sales@intactgenomics.com) | [ig@intactgenomics.com](mailto:ig@intactgenomics.com)

**Website:** [www.intactgenomics.com](http://www.intactgenomics.com)

© 2024 Intact Genomics, Inc  
All Rights Reserved

