



T4 gp32 Protein

Manual

Catalog #	3511	3512	3515	3518	3516
Package Size	100μg	200µg	500µg	1000µg	500μg
Volume	20μΙ	40μl	100μΙ	200μΙ	50μΙ
Concentration	5μg/μl				10μg/μl



Important!

-20°C Storage Required

- * Immediately inspect packages
- * Freeze upon receipt

Intact Genomics, Inc.



visit us online for more products & custom services



T4 gp32 Protein

Table of Contents

Product Description	3
Protein Purity	3
Product Source	3
Components and Storage	3
Reaction Buffer Composition	3
Heat Inactivation	3
Enzyme Concentration	4
Quality Control	5
References	5
Related Products	5
Ordering Information	5
Technical Support	6



Description:

Intact Genomics (ig®) T4 gp32 is a single-stranded DNA binding protein required for T4 DNA replication, recombination, and repair (1, 2). It improves the efficiency of reverse transcriptase (RT) during RT-PCR (3), enhances T4 DNA polymerase activity (4), as well as increases the yield of PCR products (5).

Protein Purity:

The physical purity of this enzyme is ≥98% as assessed by SDS-PAGE with Coomassie® blue staining (Fig. 1).

Product Source:

E. coli BL21 (DE3) strain expressing T4 gp32 gene.

Product Components:

- T4 gp32 protein
- 10X gp32 reaction buffer

1x GP 32 Reaction Buffer Composition:

20 mM Tris-acetate100 mM Potassium acetate 10 mM Magnesium acetate, 1 mM DTT pH 7.8 @ 25°C

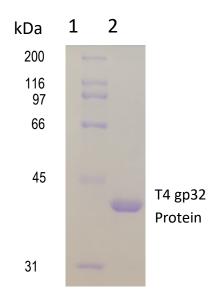


Fig. 1: Lane 1. Protein Marker Lane 2. gp32 Protein

Storage:

50 mM Tris-HCl, 50 mM KCl, 1 mM DTT, 0.1 mM EDTA 50% Glycerol, pH 7.5 @ 25°C

Heat Inactivation:

65ºC for 20 min



Enzyme Concentration:

IG uses orthogonal, 3-part approaches to determine the enzyme concentration to provide you with consistent and reliable enzymes for your needs. The quantity of a protein sample is assessed using densitometry with polyacrylamide gel electrophoresis (PAGE), UV absorbance spectra of native protein, and using a protein standard assay such as bicinchoninic acid assay (BCA) using bovine serum albumin (BSA) as a standard (Figure 1).

Why does IG use all three approaches?

- Each method above has limitations. The limitations include experimental noise, accuracy, and susceptibility to buffer and/or enzyme conditions.
- 2. Each enzyme has unique physical properties that make a single approach to analyzing proteins a challenge. Each enzyme has a different protein sequence, different requirements to be stable in solution, and different requirements to retain its maximal activity. These differences can interfere with or convolute results, especially when compared to other enzymes. When used together, however, each method provides the scientist with independent measures of both enzyme and buffer purity and quality.

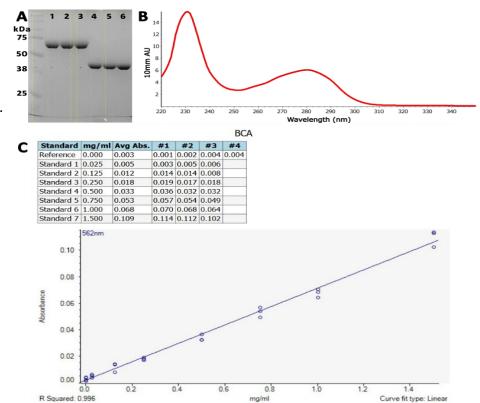


Figure 1: Enzyme quantitation methods used by IG. A) SDS-polyacrylamide gel electrophoresis. Ladder in 1st lane, 2 μ g BSA (~67 kDa) as a standard in lanes

1-3, and IG enzymes (~40 kDa) in lanes 4-6. The yellow boxes are the areas evaluated by densitometry. The integrated band intensities of IG enzymes are compared with integrated band intensities from BSA to assay concentration. B) UV spectrum of a clean IG enzyme with protein peaks at 230 nm and at 280 nm. An extinction coefficient at 280 nm is typically used to quantify protein using these spectra with buffer subtraction at 330 nm. C) BCA standard curve for BSA. The curve is used to calculate an IG enzyme concentration using BSA as the standard.



Quality Control:

T4 gp32 Protein is free from detectable nuclease activities.

References:

- 1. Chase, J. W. and Williams, K. R. (1986) Ann. Rev. Biochem. 55, 103-136
- 2. Sinha, N. K. and Snustad, D. P. (1971) J. Mol. Biol. 62, 267-271.
- **3.** Baugh, L.R. et al. (2001). Nucl. Acids Res. 29, e29.
- **4.** Topal, M.D. and Sinha, N.K. (1983). J. Biol. Chem. 258, 12274-12279.
- 5. Schwartz, K. et al. (1990). Nucl. Acids Res. 18, 1079.

Related Products:

- T4 UvsX DNA Recombinase (Cat.# 3562)
- T4 UvsY Protein (Cat.# 3572)
- Bsu DNA Polymerase (Cat.# 3585)
- Sau DNA Polymerase (Cat.# 3595)
- Exonuclease III (Cat.# 3415)
- Exonuclease IV (Nfo) (Cat.# 3425)

Ordering Information:

- Order online within the USA. Place orders on www.intactgenomics.com using our secure Shopping Cart.
- Order by email, phone, or fax.

Email: sales@intactgenomics.com

Phone: (314) 942-3655 | Toll-free: 855-835-7172 | Fax: (314) 942-3656

Order via our distributors.



Intact Genomics owns the following registered trademarks granted by the United States Patent and Trademark Office (USPTO): Intact Genomics®, IG®, ig®, igTherapeutics®, FastAmp®, i7®, DirectPlate®.

All technology protocols discussed within this manual are assumed proprietary to Intact Genomics. This Product may be covered by pending or issued patents or may have certain limitations. Please contact us for more information. Purchase of this material conveys to buyer the non-transferable right to use the material purchased in research conducted by buyer, whether for teaching, non-commercial or commercial research purposes. Buyer may not sell or otherwise transfer these materials, its components, or unmodified descendants to a third party.

Product Use Limitation and Disclaimers

This product is for research purposes only. It is not intended for therapeutic or diagnostic purposes in humans or animals. This product contains chemicals which may be harmful if misused or direct human contact is made.

Intact Genomics is dedicated to practicing and maintaining science and technology ethics. Buyer agrees to use the purchased materials in full compliance with applicable law and regulations.

Technical Support & Customer Services

Intact Genomics (IG®) is dedicated to customer satisfaction regarding the use of our products for your research needs. Each new lot of our products is thoroughly tested to ensure it meets high quality standards and provides excellent results. We appreciate your business and your feedback regarding the performance of our products in your applications. Please follow the instructions carefully and contact us if additional assistance is needed.

Our hours are Monday - Friday, 8AM to 5PM, U.S. Central Standard Time.

Intact Genomics, Inc.

11840 Westline Industrial Drive, Suite 120, St. Louis, MO. 63146, USA

Phone: (314) 942-3655 | Toll-free: 855-835-7172 | Fax: (314) 942-3656

Email: sales@intactgenomics.com | ig@intactgenomics.com

Website: www.intactgenomics.com



© 2024 Intact Genomics, Inc

All Rights Reserved